



11.6" Win10 Tablet Digital Camera, 5.0M

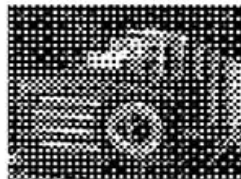



11.6" Win10 Tablet Digital Camera, 5.0M

- 11.6" LCD Resolution 1920*1080 Win10 Tablet 64 Bits
- 16M Digital Camera 1/2.3" CMOS Sony Sensor
- 2.4G/5G WIFI, Bluetooth, USB3.0, HDMI, Ethernet Port
- S-EYE Measure Software Installed
- Type C Power Supply DC12V/2A Adapter

Product Details

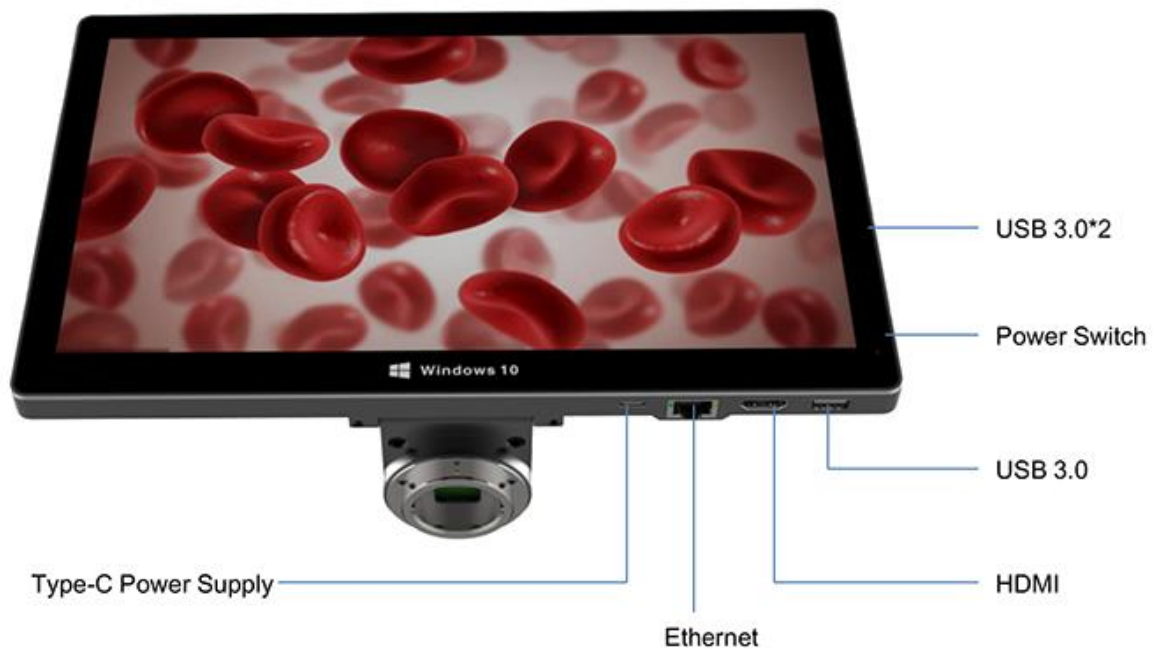

S-EYE Measure Software For Windows

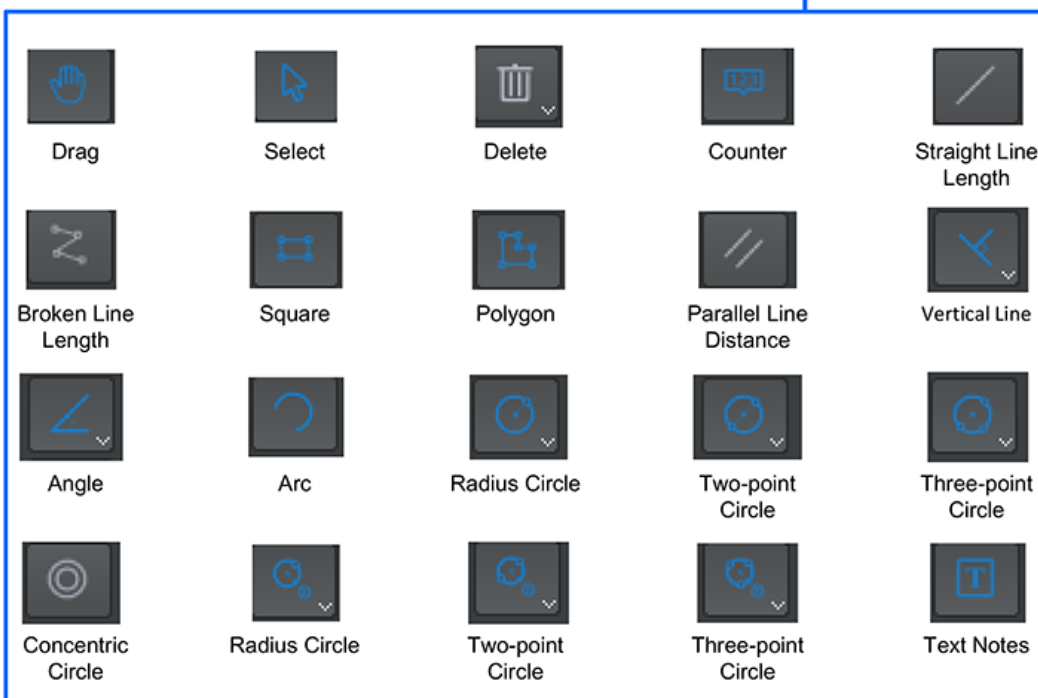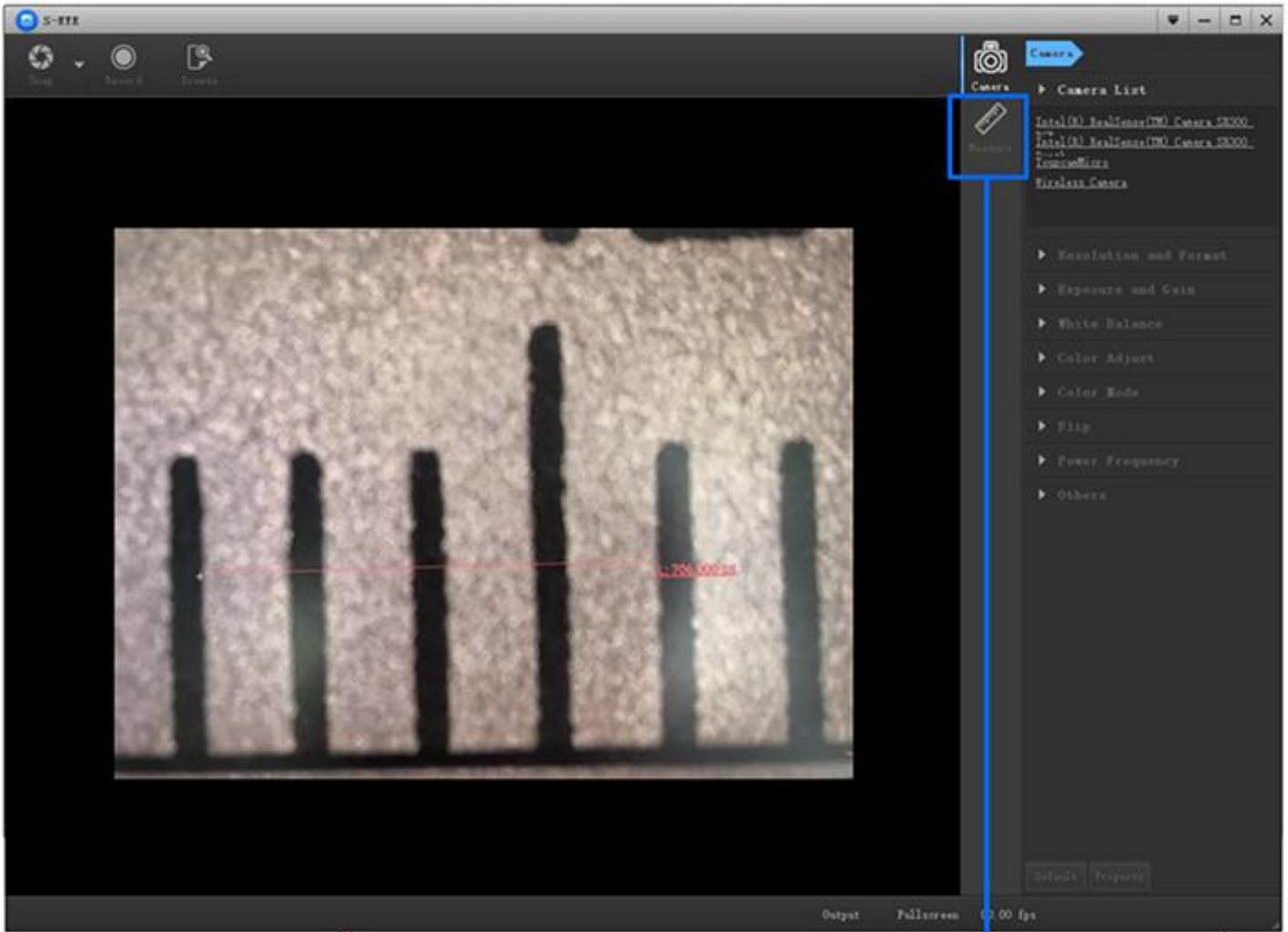



LABMICRO®

85535 45575 | 91644 42019

sales@labmicro.in / service@labmicro.in

www.labmicro.in

Technical Specification

Items	11.6" Win10 Tablet Digital Camera, 5.0M
Model No.	SLV-DT-1
Sensor	1/2.3" Sony CMOS Sensor
Resolution	16M CMOS
LCD Screen	11.6" LCD, Resolution 1920x1080p IPS Screen
CPU	Inter J4105
System	Win10 64Bits
WIFI	2.4G/5G WIFI, 802.11a/b/g/n/ac
Bluetooth	Bluetooth 4.2
Output	USB3.0, HDMI, WIFI, Ethernet, Type-C Power Supply
Pixel Size	1.335x1.335um
Power Supply	DC 12V/2A Power Adapter, Input 100-240V AC
Function	<ul style="list-style-type: none">• Photo 16M, Video 1080p Full HD• Adjust: Brightness, Contrast, Saturation, White Balance, Exposure, Flicker Suppression, Sharpness, Gamma, CB/Cr Gain• UI Interface: Mouse Control, Touch Screen• Storage Photo, Video, Preview, SD Card Format
Measure	S-EYE Measure Software: Length: Straight Line Length, Broken Line Length, Curve Length, Parallel Line Distance, Point Line Distance, Line Length Geometry: Radius Circle, Two-point Circle, Three-point Circle, Concentric Circle, Radius Circle, Two-point Circle, Three-point Circle Geometric Area: Polygon, Square Remark: Point Coordinates, Crosshairs, Coordinate Systems, Text Notes
Size / Weight	Product Size 285x177x12.5mm, Net Weight 0.9kg
Standard Package	LCD Camera, DC12V Power Adapter
Quick Manual	<ol style="list-style-type: none">1. First, connect the product to Type-C DC-12V power supply2. After connecting the power supply, press the power on button, and the blue light of the product will always be on.3. Double click the S-EYE Measure Software on the desktop of Window 104. Click "Open Video Source" in S-EYE software, and choose camera YW16006. Measure and adjust as you need.



Shop No.1, Basement floor, Best Apartments, Dr. Arunachalam Road, 5th Stage BEML Layout, Rajarajeshwari Nagar, Bengaluru - 560 098